AENOR
ER
Empresa
Registrada
ER-109/1/96

CENCOM 2

7002240

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GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.

SOMMAIRE

INFORMATION GENERALE
SOMMAIRE
LISTE D'EMBALLAGE
SPECIFICATION TECHNIQUE
DESCRIPTION DE L'EQUIPEMENT
INSTALLATION
MODE OPERATOIRE
ENTRETIEN
SCHEMA ELECTRIQUE
GARANTIE
DECLARATION DE CONFORMITE "CE"

CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
MAINTENANCE
ELECTRICAL DIAGRAM
GUARANTEE
"EC" CONFORMITY DECLARATION

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PACKING LIST

The standard equipment consist of the following components:

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Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Volume maxi. <i>Maximum volume</i>	Capacité maxi. tube <i>Maximum tube capacity</i>	Dimensions extérieures <i>Overall dimensions</i>	Vitesse maxi <i>Maximum speed</i>	R.C.F. maxi. <i>Maximum R.C.F.</i>	Régulation <i>Régulation</i>	Consom. W. <i>Consump W.</i>	Poids kg <i>Weight Kg</i>
90ml	15 ml	diam. 246 x 240 mm	3600 rpm	1780 g	8 positions <i>8 positions</i>	92	5

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EQUIPMENT DESCRIPTION

CENCOM is a benchtop centrifuge with an angle head of 6 15ml. tubes. It is supplied with safety cover and 8 speed positions.

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INSTALLATION

Place the centrifuge on a flat, horizontal, level surface.

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Do not use the apparatus to centrifuge liquids which are susceptible to form explosive mixtures.

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate of the apparatus.

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Do not use the apparatus if it is NOT EARTHED.

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If you have to change the plug bear in mind the following:

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STARTING UP:

1. Connect into mains.
2. Put the tubes into the head in a balance way. If all buckets are not used, take notice that the load must be balanced symmetrically.
3. Close the cover.
4. Select the speed. (r.p.m.):
1000, 1500, 2000, 2500, 3000, 3500, 4000.
5. Select the centrifugation time. (hh.mm)
6. Press On / Off to start centrifugation.

STOP:

Press On / Off and wait few minutes before opening the door.

Do not place tubes and remove if the head is not complete stopped.

Leave the head to stop by itself. Do not use any tool or your hand to brake the head.

SAFETY:

Microswitch which disconnects the motor when opening the cover (in case of stop omission of the motor)

CAUTION!! DO NOT OPEN THE COVER WHEN THE MOTOR IS WORKING

MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of casing, plastic and painted metals: Alcohol with cotton duster.

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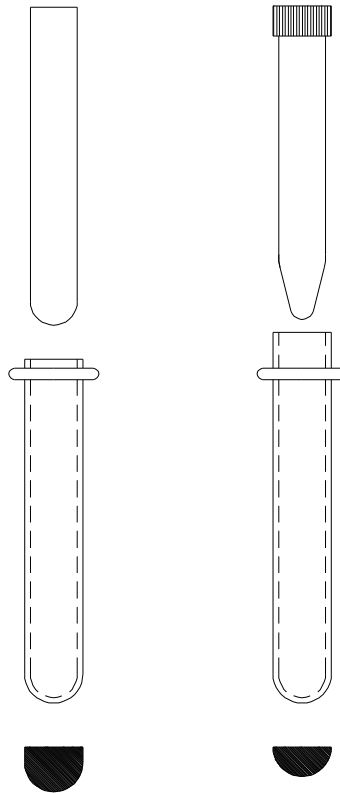
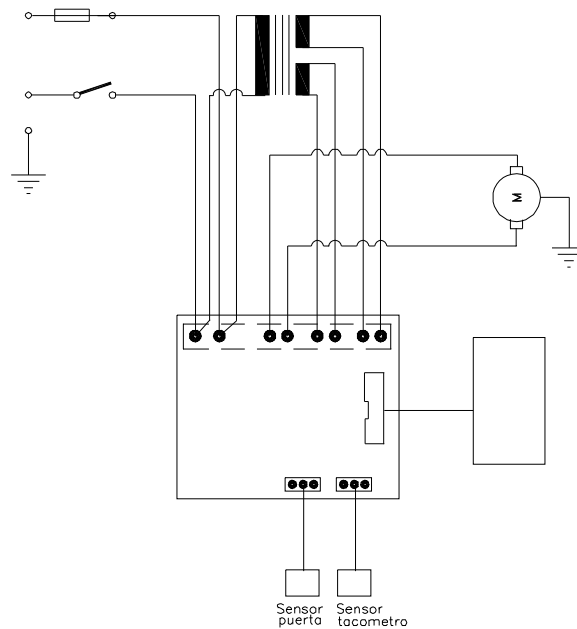
GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA,S.A..

Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA,S.A. cancels the guarantee automatically.

DECLARATION DE CONFORMITE "CE"**"EC" CONFORMITY DECLARATION**

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PROGRAMME DE FABRICATION / MANUFACTURING PROGRAMME

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