QUICK START GUIDE

EN

Quick Start Guide

Thank you for purchasing this Moticam. This camera is a delicate instrument and should be handled with care. When not in use, please store the camera in a safe place with the dust cap attached. Do not let the camera get in contact with water or expose it to extreme temperatures (below 0 or above 50 degrees Celsius). In order to receive the highest level of service, please retain your receipt as proof of purchase and **register your camera online at www.motic.com today.**

Should you have any questions please do not hesitate to **contact our customer service.**

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- I**
- <https://dl-hk.motic.com/download/moticonnect.html>

EN Using the camera on an eyepiece

1. Download the software MotiConnect:
<https://dl-hk.motic.com/download/moticonnect.html>
2. Remove the Dust Cap (B) from the camera (A) and attach the focusable lens (C).
3. Choose the right size eyepiece adapter (D) and attach this to the focusable lens.
4. Attach this assembly to your microscope's eyepiece by tightening the centering screws on the eyepiece adapter.
5. Connect the USB cable (F) onto the camera and then plug it into the computer

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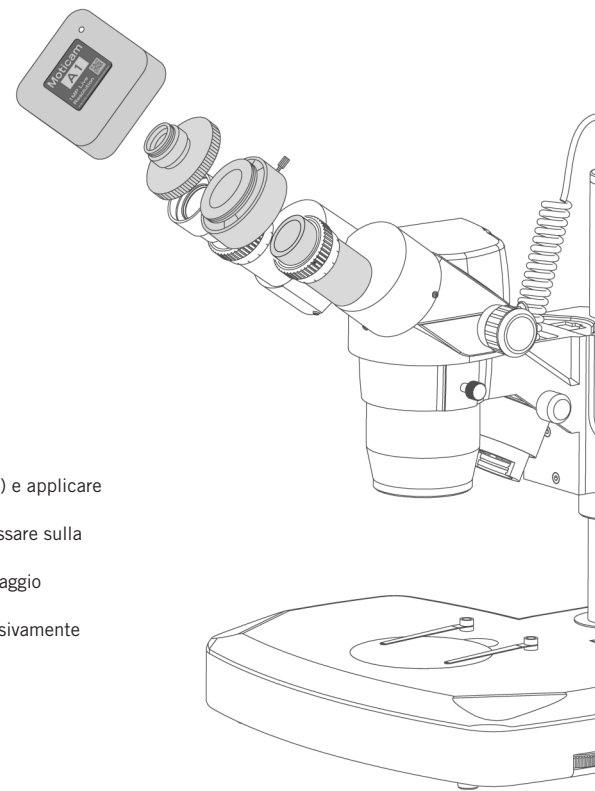
EN Using the camera as a Macro Camera

1. Download the software MotiConnect:
<https://dl-hk.motic.com/download/moticonnect.html>
2. Remove the Dust Cap (B) from the camera (A) and attach the focusable lens (C).
3. Attach the Macro Tube (H) onto the lens
4. Place this assembly over the item you wish to view and ensure there is adequate lighting. If necessary, a desk lamp is a good solution.
5. Connect the USB cable (F) onto the camera and then plug it into the computer

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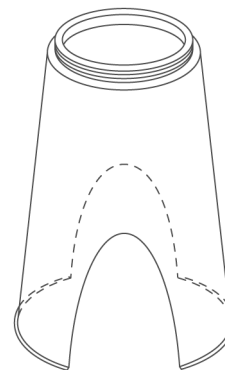
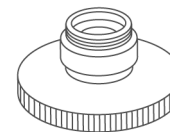
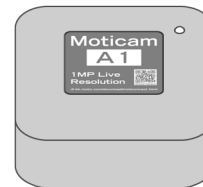
EN Using the camera on a Trinocular Microscope

1. Download the software MotiConnect:
<https://dl-hk.motic.com/download/moticonnect.html>
2. Remove the Dust Cap (B) from the camera (A) and attach the microscope's C-Mount adapter (*Please note that the C-Mount adapter is specific to your microscope. For Motic microscopes, please consult the relevant Motic Accessory list and choose an appropriate adapter for your Moticam. Depending on the type of adapter and microscope, the C-Ring (E) may be necessary between the camera and the C-mount adapter).
3. Connect the USB cable (F) onto the camera and then plug it into the computer.

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